



Mouse/Rat Testosterone ELISA

Cat. # TE187S-100 (96 Tests)

INTENDED USE

For Research Use Only. Not for use in diagnostic procedures.

MATERIALS PROVIDED	96 Tests
1. Microwell coated with Goat Anti-Rabbit IgG	12x8x1
2. Standard: 6 vials (ready to use)	0.5ml
3. Rabbit Anti-Testosterone Reagent (ready to use)	7ml
4. Assay Diluent: 1 bottle (Ready to Use)	12ml
5. Enzyme Conjugate Conc. (20X): 1 Vial	0.7ml
6. TMB Substrate: 1 bottle (ready to use)	12ml
7. Stop Solution: 1 bottle (Ready to use)	12ml
8. Wash Buffer (20X): 1 bottle	25ml

MATERIALS NOT PROVIDED

1. Distilled or deionized water. Precision pipettes. Disposable pipette tips
2. Microtiter well reader capable of reading absorbance at 450nm
3. Absorbance paper
4. Paper towel
5. Graph paper

STORAGE AND STABILITY

1. Store the kit at 2 - 8° C.
2. Keep microwells sealed in a dry bag with desiccants.
3. The reagents are stable until expiration of the kit.
4. Do not expose reagents to heat, sun, or strong light.

WARNINGS AND PRECAUTIONS

1. For Research Use Only. Not for use in diagnostic procedures.
2. For laboratory use.
3. Potential biohazardous materials:

The kit contains animal and/or human source components. All the human components have been tested and found non-reactive for hepatitis B surface antigen as well as HIV antibody with FDA licensed reagents. However, as there is no test method that can offer complete assurance that HIV, Hepatitis B virus or other infectious agents are absent, these reagents should be handled at the Biosafety Level 2, as recommended in the Centers for Disease Control/National Institutes of Health manual, "Biosafety in Microbiological and Biomedical Laboratories." 1984. All the animal products, if any, have been derived from animals of US origin and processed in USDA licensed facilities.

4. Optimal results will be obtained by strict adherence to this protocol. Accurate and precise pipetting, as well as following the exact time and temperature requirements prescribed are essential. Any deviation from this may yield invalid data.

REFERENCES

1. Chen, A., Bookstein, J.J., Meldrum, D.R., Diagnosis of a testosterone-secreting adrenal adenoma by selective venous catheterization. *Fertil. Steril.*, 1991; 55: 1202-1203.
2. Granoff, A.B. and Abraham, G.E., Peripheral and adrenal venous levels of steroids in a patient with virilizing adrenal adenoma. *Obstet. Gynecol.*, 1979; 53:111-115.
3. Bricaire, C., Raynaud, A., Benotmane, A., et al., Selective venous catheterization in the evaluation of hyperandrogenism. *J. Endocrinol Invest.*, 1991; 14: 949-956.
4. Heinonen, P.K., Androgen production by epithelial ovarian tumours in post-menopausal women. *Maturitas*, 1991; 13: 117-117-122
5. Tietz, N.W. ed., *Clinical Guide to Laboratory Tests*, 3rd Edition, W.B. Saunders, Co., Philadelphia, 1995: 578-580.
6. USA Center for Disease Control/National Institute of Health Manual, "Biosafety in Microbiological and Biomedical Laboratories"⁸⁴
7. *ICN Guide to Endocrine Testing*. Diagnostic Division, ICN Biomedicals, Inc. pp. 2:33-35; 3:4-6.

2016-06-08



- The components in this kit are intended for use as an integral unit. The components of different lots should not be mixed. It is recommended that serum samples be run in duplicate.
- Do not pipette by mouth. Do not smoke, eat, or drink in the areas in which specimens or kit reagents are handled.

SPECIMEN COLLECTION HANDLING

- Collect blood specimens and separate the serum immediately.
- Typically, specimens may be stored refrigerated at (2-8°C) for 5 days. If storage time exceeds 5 days, store frozen at (-20° C) for up to one month.
- Avoid multiple freeze-thaw cycles.
- Prior to assay, frozen sera should be completely thawed and mixed.
- Do not use grossly lipemic specimens.

REAGENT PREPARATION

- 20X Enzyme conjugate:** Prepare 1X working solution at 1:20 with assay diluent (e.g. Add 0.1ml of the Testosterone enzyme conjugate concentrate to 1.9ml of assay diluent)
- 20X Wash Buffer** Prepare 1X Wash buffer by adding the contents of the bottle (25 ml, 20X) to 475 ml of distilled or deionized water. Store at room temperature (18-26 °C).

ASSAY PROCEDURE

- Secure the desired number of coated wells in the holder.
- Dispense 25 µl of standards, specimens and controls into appropriate wells.
- Dispense 100 µl of working dilution of Testosterone-HRP Conjugate Reagent into each well.
- Dispense 50 µl of rabbit anti-Testosterone reagent to each well. Thoroughly mix for 30 seconds. It is very important to mix completely.
- Incubate at room temperature 60 minutes.
- Rinse and flick the microwells 3 times with 1x wash buffer water.
- Dispense 100 µl of TMB Reagent into each well. Gently mix for 5 seconds.
- Incubate at room temperature (20-25°C) for 15 minutes.
- Stop the reaction by adding 50µl of Stop Solution to each well.
- Gently mix 30 seconds. It is important to make sure that all the blue color changes to yellow color completely.
- Read absorbance at 450 nm with a microtiter well reader within 15 minutes.

CALCULATION OF RESULTS

- Calculate the mean absorbance value (A450) for each set of reference standards, controls and samples.
- Construct a standard curve by plotting the mean absorbance obtained for each reference standard against its concentration in ng/ml on a linear-linear graph paper, with absorbance values on the vertical or Y axis, and concentrations on the horizontal or X axis.
- Use the mean absorbance values for each specimen to determine the corresponding concentration of Testosterone in ng/ml from the standard curve.
- Any values obtained for diluted samples must be further converted by applying the appropriate dilution factor in the calculations.

EXAMPLE OF THE STANDARD CURVE

Testosterone (ng/ml)	Absorbance (450nm)
0	2.38
0.1	1.75
0.5	1.02
2.0	0.59
6.0	0.34
18.0	0.17

LIMITATIONS OF THE PROCEDURE

- Do not use sodium azide as preservative. Sodium azide inhibits HRP enzyme activities.